




# The Munro 60C

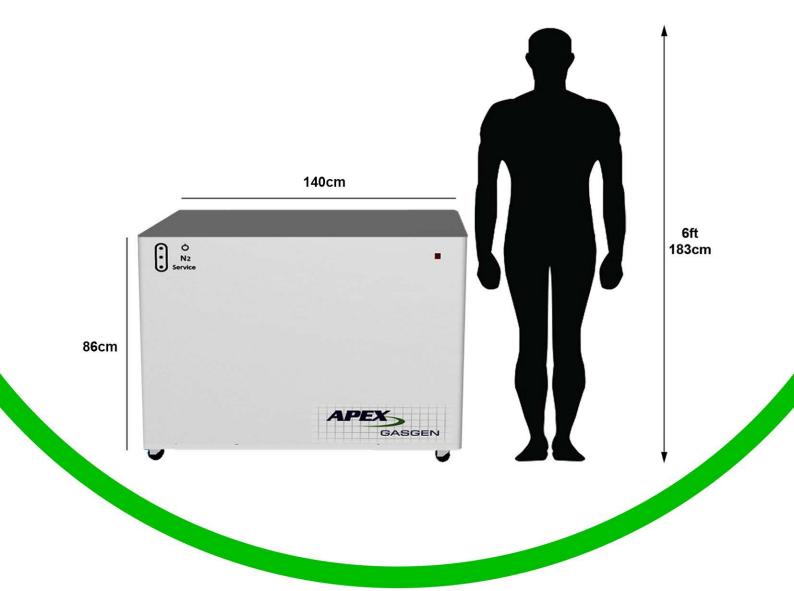

### Description

The Apex Munro 60C is a bespoke series which includes a range of models to cover most applications. When selecting a system, consideration should be given to the required flow rate, purity and delivery pressure. Additionally Apex will be happy to consult with you as to which specific model would suit you.

Featuring an integral Jun Air compressor each Nevis-C generator provides a true stand-alone solution, which offers efficency for the ideal performance time and time again.

The system uses Pressure Swing Adsorption (PSA) technology to provide you with the best possible performance. Not only are our machines highly reliable, but they are also safer to use at a better price and Apex offers to customise your system to your required specifications.

Maintenance of your Nevis Series High Purity Nitrogen Generator is simple with all regular maintenance parts being easily accessible, for peace of mind we offer a service package.

### Features & Benefits

- Standard Flow Rates up to 60 litres/min
- Output Pressure of 7 bar (100psi)
- Certified Purity Level >95%
- Medical Grade Tubing as Standard
- Stand-By Mode
- PSA Technology
- Built in Oil-Free Compressor to Reduce Contaminants
- Easy Installation
- Minimal Deterioration in Purity Over System Life Span
- Simple Maintenance and Low Cost






Model	Munro-60C
Flow / Min	60
Output Pressure (BARG/PSIG)	6.9 / 100
Dew Point (C PDP)	- 40
Power Consumption (W)	4000
Voltage (V AC)	230
Depth (mm/inch)	650 / 25.6
Height (mm/inch)	860 / 33.9
Width (mm/inch)	1400 / 55.1
Weight (kg/lbs)	200 / 441
Connection Ports	Female 1/4" BSPP
Max Altitude (m)	2000
Particulates (um)	<0.01
Lubricants	None
Humidity	80% up to 31°C decreasing linearly to 50% at 40°C

#### How It Works

Air from an external source enters the generator via a 0.01-micron filter [1] that removes contaminants and bulk moisture via an integrated auto drain. When the control system assesses that more nitrogen is required, the solenoid valve [2] is opened and pressure is allowed to build in the column [4] containing Carbon Molecular Sieve (CMS) and a residue moisture catcher [3]. This sieve adsorbs the oxygen and leaves the dry nitrogen free to flow through the non-return valve into the direction of the variable reservoir capacity (VRC) [7] Following on from this the Nitrogen is allowed to flow to the application via another 0.01-micron filter [8] and pressure and flow regulators [9]. The trapped oxygen is meanwhile filtered out through the exhaust valves [6]

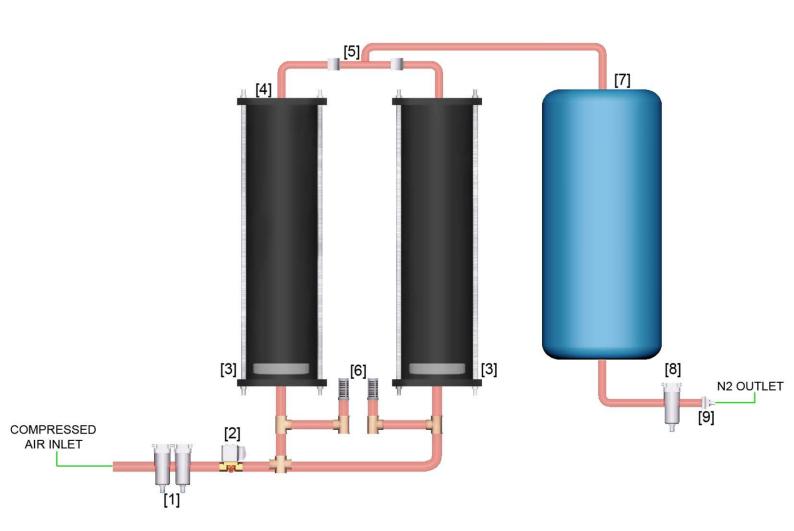



- Mass Spectrometry (LCMS instruments)

- Drying Gas & Shield Gas

- Fibre Optics

- Blanket gas for Ovens

- Special Coatings

- Wave Soldering

- Packaging

## Warranty

All Apex units are supplied with our standard 1 year warranty. Accompanied with our on-going care and servicing facilities Apex provides the highest quality peace of mind.



### **Contact Details**

Apex Gasgen Ltd
Unit 36 Elderpark Workspace
Elderpark Street
Glasgow G51 3TR
Scotland

Tel: +44 (0) 141 530 7417

Email: karen@apexgasgenerators.com

Web: www.apexgasgenerators.com

